


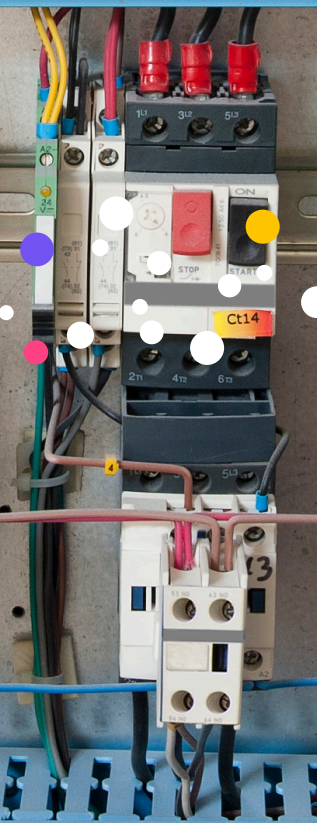

CASE STUDY :

# Wire Coatings

Results of classification of various wire coatings at different temperatures





Today, sensor networks in everything from cars to industrial manufacturing are used to ensure that the full mechanical and electrical systems are operating as intended. In industrial environments, these systems allow operators to monitor and control industrial processes to keep the full system up and running at ideal efficiency. In the case of automotive systems, the system ensures car performance and safety—sending driver alerts when a malfunction is detected.

---

In these scenarios, a heating wire can be the precursor to something as benign as a short circuit to something more catastrophic like the overheating of an entire system or assembly. Although temperature

sensors can often be used to detect extreme temperatures, being able to sense the distinct odor of a heating wire casing could enable more preventative alerts before a catastrophic failure occurs.





# Challenge

As manufacturing systems, vehicles and consumer end products rely on complex electronics systems, controls that not only regulate performance but ensure user and operator safety becomes a key focus. The challenge becomes how to integrate new sensors into an already full system—and ones that can provide meaningful, actionable data.

---

All drivers at one time or another have experienced the early warning of a smelly component or module as it warms up before it breaks. Or the very characteristic smell of an ageing brake or clutch. Many industrial operators can distinguish the smell of different components heating beyond their spec—whether it be mechanical or electrical.

Of all the existing sensors in these systems, this key parameter continues to be absent—odor sensors. Until now, e-noses have been too bulky and expensive to integrate into such an assembly. In addition, the data captured is meant for a more robust analysis—typically providing chemical intensities and breakdowns for further research. Until now, these solutions did not fit the need for a simple, compact odor diagnostic tool.





# Solution

Digital olfaction offers a better way. Digital olfaction combines biosensors, advanced optics and machine learning to mimic the human sense of smell. Digital olfaction solutions allow for simple odor detection and analysis to enable early warnings and alerts of potential malfunction.

To illustrate this, Aryballe tested two common types of wire found in these systems—18 AWG solid copper core with PVC coating and 6 AWG stranded wire with rubber insulation—to determine the ability of their NeOse Pro to classify the results.

The experiments were performed with a NeOse Pro equipped with a HeptaValve to ensure repeatability and consistency of samples. Between

two to six replicates were taken for each sample placed in 60 ml bottles, waiting two minutes between two measurements. Before measuring samples, a minimum of five minutes of waiting time was allowed to ensure complete formation of the headspace in the vial.

The experiment measurements were taken at four temperatures: ambient, 50°C, 80°C, and 150°C.

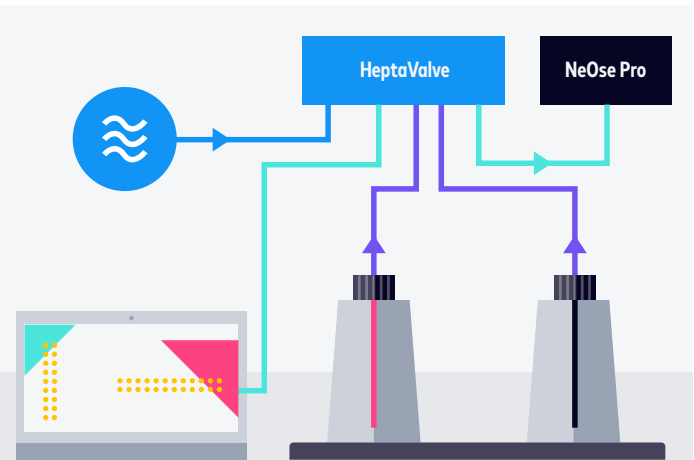


Figure 1: Experiment set up with HeptaValve and NeOse Pro.

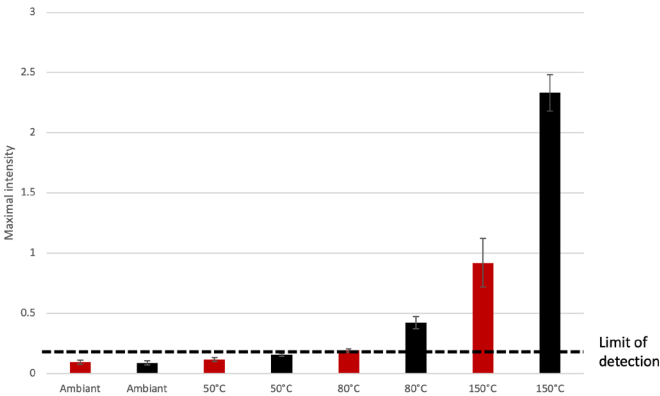


Figure 2: Intensity of headspace for two different wire samples at different temperatures.



The results of the test showed that after the five-minute waiting time, the intensity of the headspace was slightly below limit of detection for the NeOse Pro for both wires at both the ambient temperature and 50°C. Subsequent heating to 80 and 150°C provided an intensity above the NeOse Pro detection limit.

At ambient temperature, the device saw some variance in the olfactory signature replicates across samples of the same wire type (Figure 3a)—although the clustering results show two distinct zones for the wire samples (Figure 3b). Heating the samples to 50°C saw a similar variance and zone clustering, without significant improvement in the discrimination (Figure 4a and 4b).

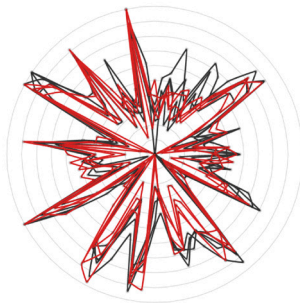


Figure 3a: Signatures of the wire samples captured by the NeOse Pro at ambient temperature.

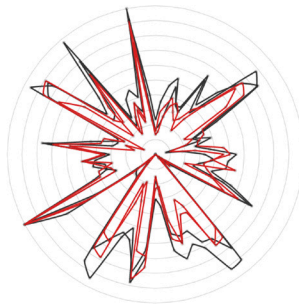


Figure 4a: Signatures of the wire samples captured by the NeOse Pro at 50°C.

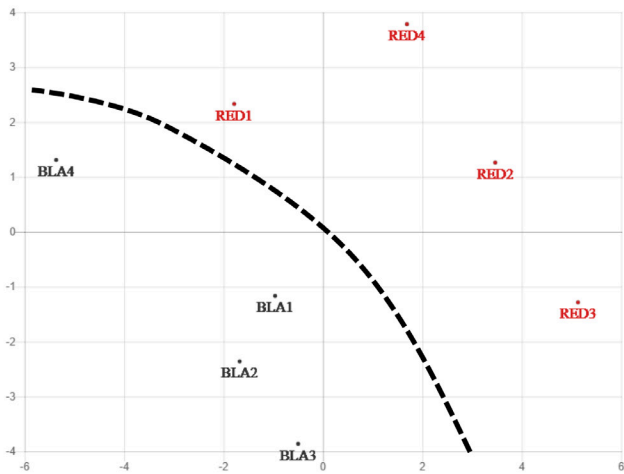


Figure 3b: PCA analysis showing zones for the red and black wires at ambient temperature.

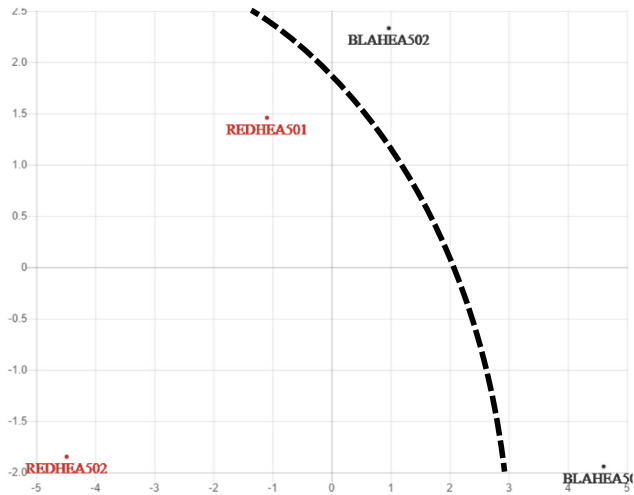


Figure 4b: PCA analysis showing zones for the red and black wires at 50°C.





As the environment continues to heat up, mimicking a heating component that begins to melt a casing, the results become more definitive. At a temperature of 80°C, there is still some variance in the olfactory signatures for the red wire, but the signatures of the black wire start to become more uniform (Figure 5a). Since the intensity of the wire at 80°C is over the NeOse detection limit, we start to see two zones with significant distinction in the clustering results (Figure 5b).

Finally, when the temperature was increased to 150°C, we see reproducible olfactory signatures across all replicates for both the red and black wires (Figure 6a) with a significant distinction in the mean olfactory signatures across the products (Figure 6b).

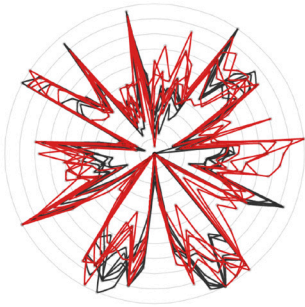


Figure 5a: Signatures of the wire samples captured by the NeOse Pro at 80°C.



Figure 6a: Signatures of the wire samples captured by the NeOse Pro at 150°C.

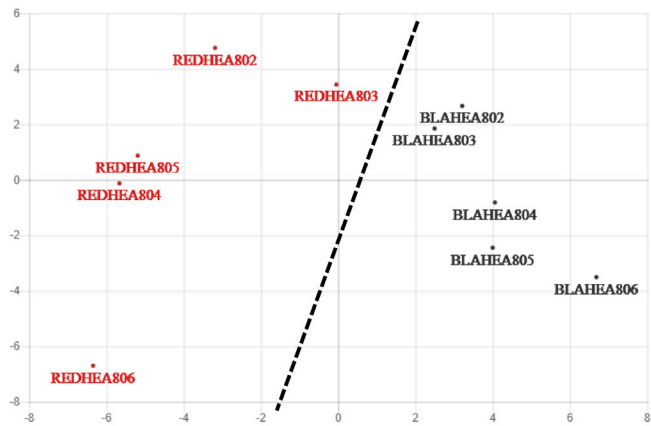


Figure 5b: PCA analysis showing zones for the red and black wires at 80°C.

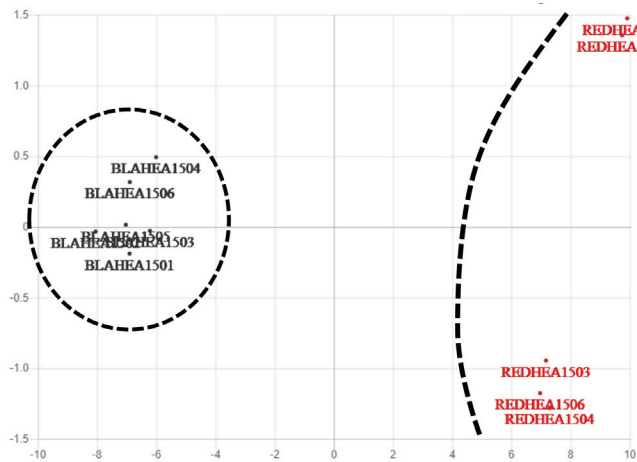


Figure 6b: PCA analysis showing clustering for the red and black wires at 150°C.





# Conclusion

The results of the tests showed that Aryballe's digital olfaction technology can distinguish between wire types at high temperatures—those thought to signify component overheating or system failures. Additional studies between temperatures from 80 to 150°C could be performed to determine the exact temperature where a consistent discrimination can first be seen.







## About Aryballe

The best human noses can distinguish 10,000 odors. Unlike color and sound, smell does not fall along a clear spectrum, making it hard to compare various odors. But Aryballe is helping to change this by evaluating characteristics of individual scent molecules and testing them against a data-base of known smells using a combination of biochemistry, advance optics and machine learning.

Based in Grenoble, France, Aryballe develops and manufactures bio-inspired "digital nose" sensors enabling groundbreaking applications in the food, cosmetics and automotive industries. Founded in 2014, it released its first product, the digital nose NeOse Pro in early 2018. Fast, portable and sensitive to hundreds of odors, NeOse Pro is used for quality control in the cosmetic industry, new flavors development in the food & beverage industry, or materials quality monitoring in the automotive industry.

---

### HEADQUARTERS

7 rue des Arts et Métiers  
38000 Grenoble, France  
+33 4 28 70 69 00

### US

101 Crawfords Corner Rd  
Suite 4-101R  
Holmdel, NJ 07733 USA

 [aryballe.com](http://aryballe.com)



© 2019 Aryballe Technologies. All rights reserved. Aryballe, NeOse, and the Aryballe logo are trademarks of Aryballe Technologies and/or its affiliates. All other trademarks are the property of their respective holders.